



MMM Beverage
Partner Kft.

Detection and identification of viable spoilage Yeasts and Moulds

FUN LC RT-PCR Workflow System

MMM Beverage Partner's FUN LC RT-PCR Workflow System allows rapid detection of viable spoilage yeasts and moulds from filterable and non-filterable beverage products (juices, teas, energy drinks, concentrates). Relevant yeasts and moulds can be distinguished e.g. *Candida parapsilosis*, *Candida pseudointermedia*, *Debaryomyces hansenii*, *Dekkera spp.*, *Metschnikowia pulcherrima*, *Saccharomyces cerevisiae*, *Zygosaccharomyces spp.*, *Yarrowia lipolitica*, *Aspergillus spp.*, *Mucor spp.*, *Paecylomyces spp.*, *Penicillium spp.*, *Rhizopus spp.*, *Talaromyces spp.*, *Trichoderma spp.*, *Ulocladium spp.* (Fig 1.). Detection and identification is carried out in one step.

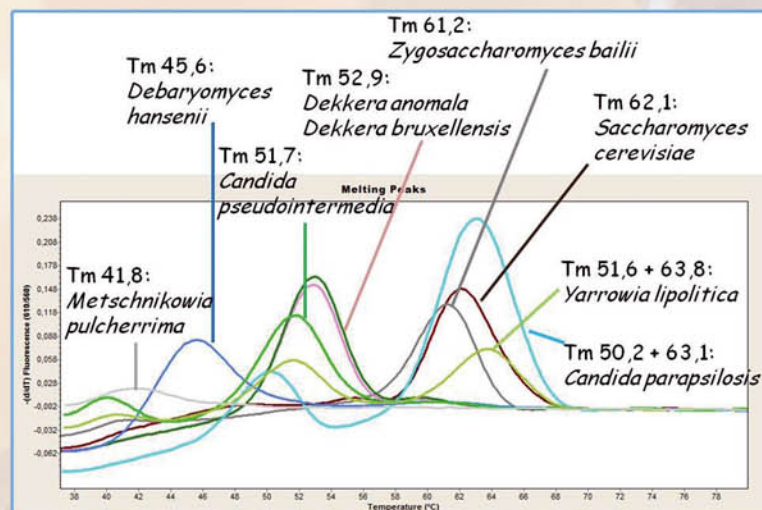


Fig 1. Differentiation of yeasts

MMM Beverage Partner's FUN LC RT-PCR Workflow System uses automated nucleic acid purification, automated PCR set up and automated detection and identification in results reporting.

The FUN LC RT-PCR Workflow System Kit includes:

- FUN Pre-cultivation Kit
- FUN Isolation Kit
- Yeast and Mould detection Real-Time PCR Kit.

FUN Pre-cultivation Kit is specifically developed for yeast & mould cultivation, enzymatic treatment and cell-disruption for non-filterable and filterable samples prior to nucleic acid Isolation in FUN LC RT-PCR Workflow System. Yeasts & moulds are cultivated at selective conditions. After the pre-cultivation the free nucleic acids are eliminated then the intact cells are disrupted prior to nucleic acid isolation.

FUN Isolation Kit is designed for the purification of genomic DNA of fungi from non-filterable and filterable samples using the MagNA Pure LC Instrument (Roche). The purified high molecular weight DNA is used for yeast & mould detection Real-Time PCR Kit.

Microbiological Quality Control of Beverages

Rapid Detection and Identification of Viable Spoilage

Microorganisms

Yeast and Mould detection Real-Time PCR Kit is specifically developed for real-time, online PCR using Hybridization Probes. Yeast & mould DNA is amplified with specific primers designed for the conservative region of the fungi genome. The amplicon is detected by fluorescence using a specific pair of Hybridization Probes. Hybridization Probes are used for distinguishing of the different yeast & mould species.

Sample material:

There is no need to separate any sample components (e.g. fruit fibers) before the analysis which separation would mean the risk to lose contaminants attached to the separated components. Our system provides you the possibility to analyze the total contents of your sample even if it is 100% fruit juices with pulp.

- bottled final beverage products (filterable and non-filterable, including 100% fruit juices with pulp)
- concentrates
- pure culture (e.g., colonies)
- water samples
- swab samples

Key features of our system:

- Fast: Result can be gained even within 3 hours.
- Robust: 100ml can be examined even from non-filterable beverages.
- Throughput of the LC RT-PCR Workflow System: 60 bottles / 1.5 h
- Detection and identification of the contaminants in one step
- Sensitive: Even the presence of one single cell of unwanted microorganism in the sample can be detected and identified.
- Specific: While all of spoilage microorganisms are detected, harmless microorganisms out of interest do not affect the result due to specific hybridization probes.
- The same kit and protocol can be used for filterable and for non-filterable samples.
- Automated processes
- Automatic evaluation and interpretation of results
- Electronic storage of data
- Effective tool for both Quality assurance and Root Cause analysis
- Flexible system
- Reproducible results

We are proud of our Real-Time PCR system being the first rapid PCR based microbiological method approved and recognized by Coca-Cola Hellenic as an alternative method to the conventional analysis. Coca-Cola Hellenic introduced our Real-Time PCR systems in their quality assurance system for detection of potential spoilage microorganisms.

Related products:

- **ACB LC RT-PCR Workflow System Kit:** Rapid detection of viable thermo-, acidophilic bacteria belonging to the genus of *Alicyclobacillus* and identification of different *Alicyclobacillus* species from filterable and non-filterable beverage products (juices, teas, energy drinks, concentrates).
- **LAB LC RT-PCR Workflow System Kit:** Rapid detection of viable *Lactobacillus* bacteria from filterable and non-filterable beverage products (juices, teas, energy drinks, concentrates). Coming soon...
- **RT-PCR upgrade package:** This product allows you to upgrade your classical microbiological laboratory to RT-PCR laboratory. This upgrade package includes all the tools you need: equipments, specific software parts for nucleic acid purification and detection, installation and training.

Contact information...

Mailing address

MMM Beverage Partner Ltd.
Otthon u. 28/b.
1118 Budapest
Hungary

Laboratory

MMM Beverage Partner Ltd.
Építők útja 2-4.
2040 Budaörs
Hungary
Phone: +36 23 428896
E-mail: info@mmmbeverage.hu
web: www.mmmbeverage.hu